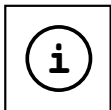
- DE Bedienungsanleitung**
- EN Operating Instructions**
- FR Mode d'emploi**
- NL Handleiding**
- IT Istruzioni per l'uso**
- ES Instrucciones de uso**
- PT Manual de utilização**

MANUAL DOWNLOAD:

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MICROSCOPE GUIDE:

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MICROSCOPE FAQ:

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EXPERIMENTS:

www.bresser.de/downloads

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- **Choking hazard** – This product contains small parts that could be swallowed by children. This poses a choking hazard.
- **Risk of electric shock** – This device contains electronic components that operate via a power source (power supply and/or batteries). Only use the device as described in the manual, otherwise you run the risk of an electric shock.
- **Risk of fire/explosion** – Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- **Risk of chemical burn** – Make sure you insert the batteries correctly. Empty or damaged batteries could cause burns if they come into contact with the skin. If necessary, wear adequate gloves for protection.
- Leaking battery acid can lead to che-

mical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.

- Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. Remove the batteries from the unit if it has not been used for a long time.
- Never recharge normal, non-rechargeable batteries. This could lead to explosion during the charging process.
- Rechargeable batteries may only be charged under adult supervision.
- Rechargeable batteries must be removed from the toy before charging.
- Do not short-circuit the terminals.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Ser-

vice Centre and can send the device in to be repaired, if necessary.

- Tools with sharp edges are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- Keep instructions and packaging as they contain important information.

DISPOSAL

Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2012/19/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

Empty old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 01.06.2006 from your local waste disposal service or environmental authority.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please pay attention to dispose of your used batteries as required by law - at a local collection point or in the retail market (a disposal in domestic waste violates the Battery Directive).

Batteries that contain toxins are marked with a sign and a chemical symbol. „Cd“ = cadmium, „Hg“ = mercury, „Pb“ = lead.

Cd¹

Hg²

Pb³

¹ battery contains cadmium

² battery contains mercury

³ battery contains lead

EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address: www.bresser.de/download/8852001/CE/8852001_CE.pdf

UKCA Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the UKCA declaration of conformity is available at the following internet address: www.bresser.de/download/8852001/UKCA/8852001_UKCA.pdf

Bresser UK Ltd. • Suite 3G, Eden House, Enterprise Way, Edenbridge, Kent TN8 6HF, Great Britain

Your microscope is made up of the following parts (Fig. 1-4):

- 1 Binocular view piece (with eyepiece supports)
- 2 Stereo objective
- 3 Eyepieces
- 4 Illumination (LED)
- 5 Microscope stage (with object plate)
- 6 Stage clips
- 7 Focus knob
- 8 On/off switch
- 9 Battery compartment
- 10 Mounting screws
- 11 Battery contacts (flat sheets and spring terminals)

12 mineral samples

Incident light microscopy

5 prepared slides

Transmitted light microscopy

Object plate

Location

Before you start, choose an ideal location for using your microscope. It is important that you have enough light available (window, desk lamp).

Simple Observation

Place the object to be observed (for example, a rock) in the middle of the microscope stage (5). When the object is large enough, it is best to fix it in place with the stage clips (6).

Now, take a look through the binocular eye lens (1) and change the distance of the eyepiece (3), so that you have a circular, clear image.

The sharpness is adjusted by slowly turning the focus knob (7).

Illumination

Remove both eyepieces (3) from the eyepiece supports (1) and turn the entire microscope over. In the base of the microscope there is a battery compartment (9). Loosen the fixing screw (10) of the battery compartment cover and remove the cover.

Place the batteries (2x AA) in the compartment so that the flat minus poles (-) press against the spring terminal and the plus poles (+) are touching the flat contact sheets (11).

Put on the cover and fix it with the screw (10). Turn the microscope over again. Place the eyepieces back on. You can not turn on the light (4) with the on/off switch (8) on the base of the microscope.

Experiments

Use the following web link to find interesting experiments you can try out.

How do you hatch brine shrimp? How to make salt crystals and much more.

<http://www.bresser.de/downloads>

TIPS on cleaning

Remove the device from its energy source before cleaning (remove batteries).

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

Clean the lens (objective and eyepiece) only with a soft lint-free cleaning cloth (e.g. micro-fibre). Do not use excessive pressure - this may scratch the lens.

Dampen the cleaning cloth with a spectacle cleaning fluid and use it on very dirty lenses.

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